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於 N. M. I. M. R.

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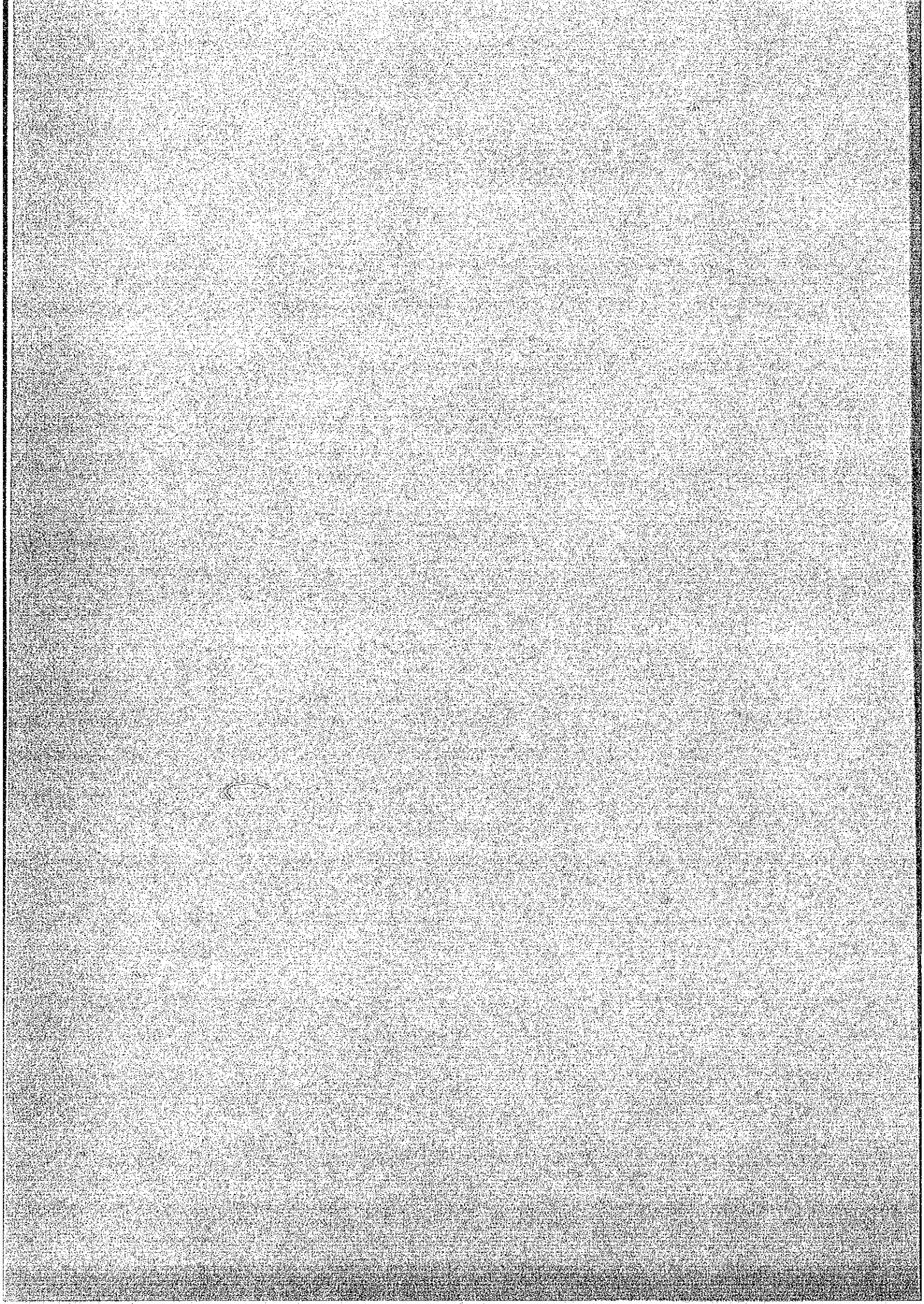
N.M.I.M.R

Republic of Ghana

Re.: Result of Inspection and Repair of Equipment for the
Noguchi Memorial Institute for Medical Research (N.M.I.M.R.)

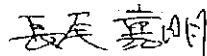
Dear Nkruma,

We wish to inform you that with your workshop staff's cooperation, our inspection and repairing work has been successfully completed. Therefore, we are submitting the following Report on Repair of your Equipment. Please find attached the details of our report.

We very much appreciated your kind cooperation during our stay in Ghana.

Thanking you, we are,

Sincerely yours,

Yoshiakira NAGAO

Leader

Medisun Corporation

1. Member of Repair Team

- | | | |
|----------------------|----------|---------------------|
| (1) Yoshiakira NAGAO | : Leader | Medisun Corporation |
| (2) Katsunori ISOBE | : Member | Medisun Corporation |
| (3) Tamotsu NOZAKI | : Member | Medisun Corporation |

2. Period of Work in Ghana

- (1) First Stage : September 20, 1993 to September 25, 1993 (6 days)
(2) Second Stage : January 10, 1994 to January 20, 1994 (11 days)

3. Scope of Work

- (1) Adjustment and repair of equipment which as requested from N.M.I.M.R.
(Refer to the Annex-1 : Summary of Inspection and Repair Report)
- (2) Inspection of the existing equipment in N.M.M.I.R.
- (3) Research of the local agents (for Generator, Air conditioner, ...etc.)
- (4) Instruction in repairing and maintenance technics to the N.M.M.I.R. staff
- (5) Workshop Programme (maintenance lectures)
(Refer to the Annex-2 : Summary of Workshop Programme)

Result of Inspection/Repair

No	Equipment	Manufacturer	Model	Condition of Equipment	Procured Parts	Qty	Result of Inspection/Repair
A1	High Speed Centrifuge	Hitachi	80P-7	Burnt out drive unit, leaking vacuum pump	Nil		Adjusted (completely)
A2	Electron Microscope	Hitachi	H-600	Defective CPU board, replace defective parts	1) Ampere meter 2) Meter, for PN-1 3) Filament 777-0179 4) O-ring L458844 5) Pressure hose 6) Pressure hose 7) Oil MR100 8) Freeze oil 56SII 9) Oil mist trap (L) 10) Fluorescent screen 11) Fluorescent screen	1 1 1 3 4 4 10 10 2 1 1	Repaired (completely)
1	Amino Acid Analyzer	Iriwa Niki	A-3300	The reactor does not work, it attains the set temp., but does not start the buffer pump	Nil		Repaired (completely)
2	Electronic Balance	A & D	EV-60A	Defective CPU board, unrepairable, replace by new one	Procured new one	1	(A & D Model: FY-21)
3	Ice Maker	Ebara	RF-51-SC	Good condition	Nil		Adjusted (completely on 1st stage)

COUNTRY Ghana
PROJECT N.M.M.I.R.

No. 2/7

Result of Inspection/Repair

No	Equipment	Manufacturer	Model	Condition of Equipment	Procured Parts	Qty	Result of Inspection/Repair
4	pH Meter	Horiba	S769-9	Damaged electrode, replace with new one	1) Electrode 5329-10C	1	Repaired (completely)
5	Deep Freezer	Ebara	ESL-260A	Faulty compressor, unrepairable	Nil		Dispose
6	Safety Cabinet	Dalton	BSC-1300	Suction device was not functioning	1) UV lamp (15W) 2) Glow lamp 3) Foot switch 4) Pipet aid PA-200	10 10 2 1	Repaired (completely)
7	Drying Oven	Sakura	TK-21	Thermo regulator sensor was defective	Nil		Repaired (completely on 1st stage)
8	Microscope	Olympus	B12	Phase contrast attachment needs to be replaced	1) Phase contrast attachment B12-PC	1	Repaired (completely)
9	Atomic Absorption Spectrophotometer	Shimadzu	UV-630-2	Defective arsenic cathode lamp	1) Arsenic cathode lamp	2	Repaired (completely)
10	Centrifuge	Sakura	300S-1	Faulty speed control	Nil		Adjusted (completely on 1st stage)
11	Spectrophotometer	Shimadzu	UV-120-02	Digital panel meter fluctuates	1) Quartz sel 10ma 2) UV lamp 3) Tungsten lamp	4 2 2	Repaired (completely)
12	Deep Freezer	Rebuko	UL11635-CH	Burnt compressor, unrepairable			Dispose

Result of Inspection/Repair

No	Equipment	Manufacturer	Model	Condition of Troubles	Procured Parts	Qty	Result of Inspection/Repair
13	Refrigerator	Hitachi	R-4867D	Defective condensing unit, unrepairable	Nil		Dispose
14	Refrigerator	Sanyo	SR-480F	Burnt compressor, unrepairable	Nil		Dispose
15	Deep Freezer	Rebuko	ULT-1285-CFD	Burnt compressor, unrepairable	Nil		Dispose
16	Refrigerator	Hitachi	R-252H	Burnt evaporative fan motor, unrepairable	Nil		Dispose
17	Refrigerator	USA GE	CAF160B	Burnt compressor, unrepairable	Nil		Dispose
18	Refrigerator	USA GE	CAF160VC	Burnt compressor, unrepairable	Nil		Dispose
19	Refrigerator	Ebara	UF220	Burnt compressor, unrepairable	Nil		Dispose
20	Refrigerator	USA GE	CAF160VC	Burnt compressor, unrepairable	Nil		Dispose
21	Freezer	Sanyo	SCR-350	Burnt compressor & condensor fan motor	Nil		Dispose
22	Refrigerator	Hitachi	RS-2603	Burnt compressor, unrepairable	Nil		Dispose
23	Steam Sterilizer	Sakura	FRA-G09	Defective sequence of operation			Repaired (completely on 1st stage)
24	Deep Freezer	Ebara	UF-22	Burnt compressor, unrepairable	Nil		Dispose

Result of Inspection/Repair

COUNTRY Ghana

PROJECT N.M.M. I. R.

No. 4/7

No	Equipment	Manufacturer	Model	Condition of Equipment	Procured Parts	Qty	Result of Inspection/Repair
A-1	Microscope	Olympus	WANOX	Defective lamp & lamp holder	1) Lamp 12V100W HAL-L 2) Lamp holder A2-LS	10 1	Repaired (completely)
A-2	Air Conditioner	Sanyo	SPR-L200	Faulty compressor (5 units)	1) Compressor	5	Recommend local procurement
A-3	Parts for Refrigerator				1) Extension valve 2) Thermostat	6 3	(BHX-56140BH...3, others...3)
A-4	Parts for Repair Tools				1) Vacuum pump 2) Vender (5/8", 3m/m)	1 2	(each 1 pc.)
A-5	Generator	Medensha	ZK-187LB	Unrepairable	Nil		Recommend local procurement
A-6	Autoclave	Sakura	ASV-3001	Defective timer	1) Timer 2) Exhaust valve knob	1 1	Repaired (completely)
A-7	Autoclave	Sakura	AC-370AE	Defective pressure control device	1) Pressure control 2) Door knob	1 1	Repaired (completely)
A-8	Parts for Balance	A & D		Adaptor for 12V	1) Adaptor (100-12V)	1	
A-9	Parts for Incubator	MPCO		Defective CO ₂ regulator (for 5 units)	1) CO ₂ regulator	5	
A-10	Parts for Water Still	Yamato	WA-550		1) Ion exchange resin	1	

COUNTRY Ghana
PROJECT N.M.M.I.R.

No. 5/7

No	Equipment	Manufacturer	Model	Condition of Equipment	Procured Parts	Qty	Result of Inspection/Repair
A-11	Light Source for Microscope	Olympus	LSK	Defective light source	1) Light source (240V 20WGE)	2	Repaired (completely)
A-12	Chromatograph Recorder	Shimadzu	C-867A	Defective recording roller			
A-13	Weighing Scale (100kg)	Yamato		Defective glass	1) Glass 2) Flame for glass	2	Repaired (completely)
A-14	Health Meter (100kg)	Yamato		Defective main body, unrepairable replace new one	Procured new one	1	
A-15	Grinder	Hitachi	GR-21	Defective switch box	1) Switch box	1	Repaired (completely)
A-16	Parts for Work Shop				1) Gas cylinder	1	Empty cylinder, for 6kg
A-17	Water Still	Yamato	WGA-28	Good condition, instruction in daily maintenance	Nil		Adjusted (completely on 1st stage)
A-18	Plate Reader	Nihon Saiyaku	Uniskan 1	Request instruction manual	1) Instruction manual	1	
A-19	Ultrasonic Homogenizer	Tomy	UR-200P	Request instruction manual	1) Instruction manual	1	
A-20	Deep Freezer	Revco	UL-T-1285C	Burnt condenser, unrepairable			Dispose

COUNTRY Ghana
PROJECT N.M.M.I.R.

No. 7/7

Result of Inspection/Repair
(Request parts from Work Shop in N.M.I.R.)

No	Equipment	Manufacturer	Model	Condition of Equipment	Procured Parts	Qty	Result of Inspection/Repair
1	Refrigerator	Amesha	P60345-724		1) Compressor 2) Dryer 3) Freon 12 4) Freon 22 (kg) 5) Copper coil 5φ, 8φ	1 2 50 10 4	Impossible procurement in Japan " " " each 2
2	Clamp for Tester				1) Clamp	1	
3	Fuse				1) 0.5, 1.0, 3.5 A	30	each 10
4	Grinding Stone	Hitachi	GR21		1) Grinding stone	10	
5	3 Wire Extension Cable				1) 1.6 x 100 m	2	(with ground)
6	2 Wire Extension Cable				1) 1.6 x 100 m	2	(Standard)
7	Vehicle	Suzuki	SJ50		1) Handle door RH 2) Handle door LH 3) Filter fuel 4) Pump assembly fuel	2 2 4 2	
8	Vehicle	Toyota	BB21		1) Filter fuel 2) Assembly fuel filter 3) Cylinder assy. (RWB)	4 2 4	

Workshop Programme (Maintenance Lectures)

Date : January 18, 1994

Objective Equipment : High Pressure Steam Sterilizer (Sakura)

Contents : How to trouble-shoot and temporarily repair
How to carry out routine maintenance and management

Attendants : N.M.M.I.R.
Mr. S.K.A. Jones
Mr. J.A. Kortei
Mr. Asiedu-Acheampong
Mr. E.O. Lamptey
(others-2 engineers)

Summary of Lectures

(1) How to maintain equipment without breakdowns for a maximum amount of time.
The operator should thoroughly understand the equipment and observe the following;

- 1) The purpose of the equipment
- 2) How to handle and operate
- 3) The instruction manual and operation manual should be well read before use
- 4) Always clean the equipment before and after use
- 5) Always check the condition of the equipment
- 6) Keep the area where the equipment is installed clean at all times
- 7) Check and control the environmental conditions (Temperature, humidity,

..., etc.) in the room where the equipment is installed

If the above mentioned conditions are strictly observed, the equipment will work effectively for a maximum amount of time.

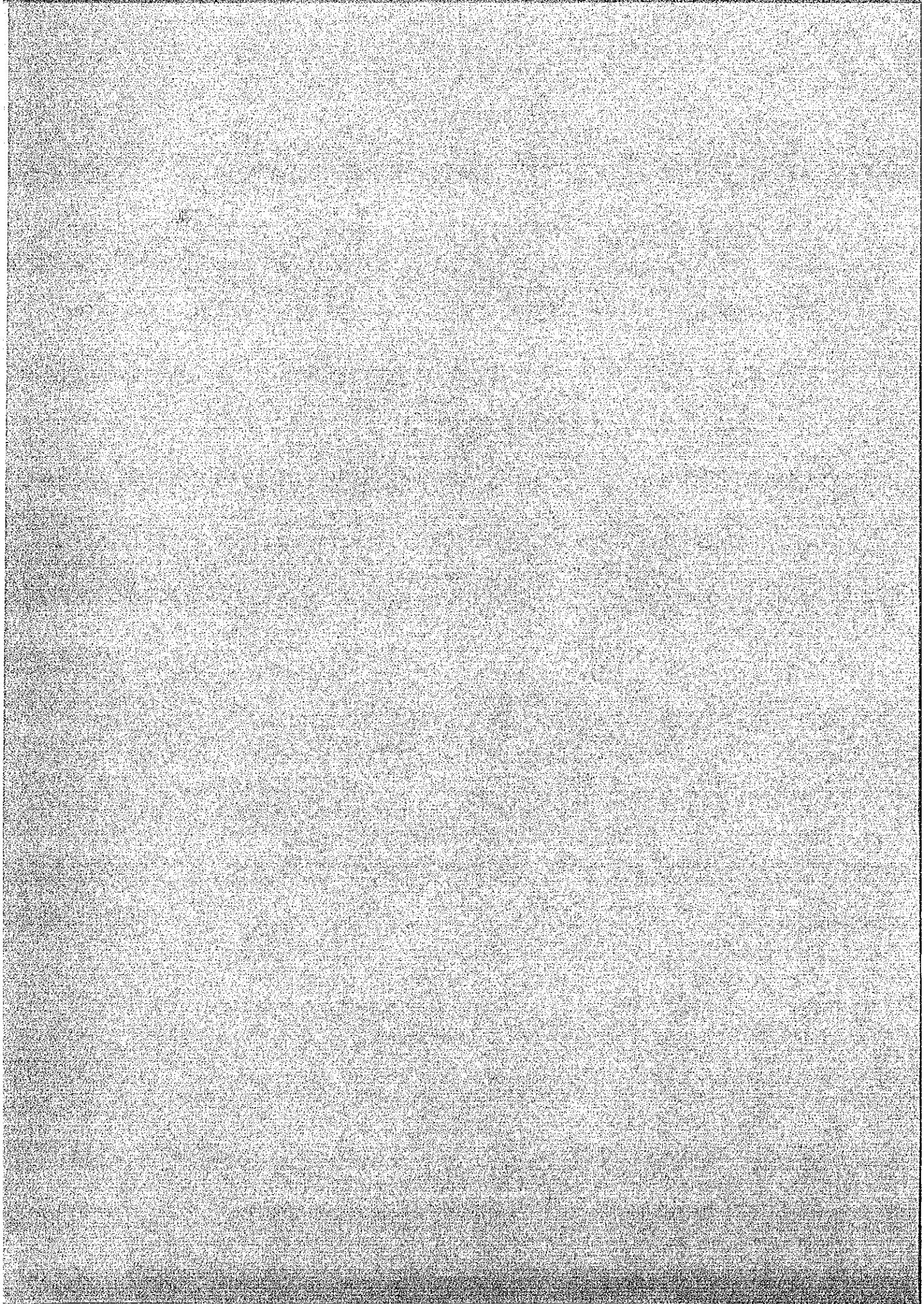
(2) The maintenance engineers should have technical knowledge of the equipment; for example the construction and how to operate, operation theory, replacement of parts and preventive maintenance.

The maintenance team should observe the following:

- 1) The engineers should always show the operator how to handle the equipment properly
- 2) Establish a routine maintenance schedule and observe that it is carried out as stipulated on maintenance chart

We hope that the maintenance engineers will be able to more effectively carry out their work after observing the above mentioned points.

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RS.8/247.

NOGUCHI MEMORIAL INSTITUTE
FOR MEDICAL RESEARCH
P. O. BOX 25
LEGON

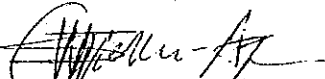
21st September, 93

Dear Sir/Madam,

JICA REPAIR TEAM - 1993

Attached is a record of the discussions that took place between the Institute and the above team on 20th September, 1993, for your information.

Yours faithfully,

S.W. Opoku-Agyakwa
ADMINISTRATIVE SECRETARY

Distribution

The Director
Participants.

MEETING WITH THE JICA REPAIR TEAM ON 20TH
SEPTEMBER, 1993 AT THE CONFERENCE ROOM

It is a policy under the Ghana-Japan Medical Co-operation Agreement that every three years, JICA sends a repair team to inspect plant and equipment donated under the agreement to evaluate their state of repair and performance, and recommend remedial action.

Such a repair team arrived in the country on 19th September 1993 to carry out the triennial exercise.

The Institute met the team on 20th September at the Conference Room to exchange ideas.

The following emerged from the meeting:

- i) the mission of the team was to take census of spare-parts required for the equipment that had broken down and submit the necessary report to the JICA Headquarters. The team would return to the country to carry out actual repairs in January 1994;
- ii) previous reports of the Institute would be the point of reference of the inspection. Thereafter the team would examine and consider requests from staff with the view to determining repair and spare-parts needs of other facilities outside the list in hand;
- iii) a copy of the report of the team at the end of the exercise would be given to the Institute.

The following were present at the meeting:

Meeting with the JICA Repair Team

1. Dr. E.A. Afari - Epidemiology Unit (in the Chair)
2. Mr. Tamotsu Nozaki - Repair Team Member
3. Mr. Yoshiakira Nagao - Leader, Repair Team
4. Mr. Katsonori Isobe - Repair Team
5. Dr. K. Taniguchi - JICA Team Leader
6. Dr. Takuro Arishima - JICA Expert
7. Dr. Y. Wagatsuma - JICA Expert
8. Dr. T. Rikimaru - JICA Expert
9. Dr. Y. Kamiya - JICA Expert
10. Mr. A.K. Ayim - E.M. Unit
11. Mr. L. Asiedu-Acheampong - Maintenance Unit
12. Mr. Jonas Asigbee - Parasitology Unit
13. Mr. Edward Addo - Nutrition Unit
14. Mr. E.O. Lamptey - Maintenance Unit
15. Mr. J.A. Kortei - Maintenance Unit
16. Mr. S.K.A. Jones - Maintenance Unit
17. Ms. J. Yartey - Nutrition Unit
18. Mr. M.A. Appawu - Epidemiology Unit
19. Dr. M. Osei-Kwasi - Virology Unit
20. Dr. N-A. Ankrah - Clinical Pathology Unit
21. Dr. M.E. Aryeetey - Parasitology Unit
22. Dr. Patience Akpedonu - Bacteriology Unit
23. Mr. S.B. Owusu - Bacteriology Unit
24. Mr. S.W. Opoku-Agyakwa - Secretary.